Microscopio Metalográfico, modelo 101 | Metallurgical microscope, model 101

Marca|Trademark:Zuzi

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Microscopio Metalográfico, modelo 101 | Metallurgical microscope, model 101

Marca|Trademark:Zuzi

Zuzi metallurgical microscope model 101 has been especially conceived for observation of any kind of opaque surfaces such as metals, alloys or mineral, by means of the incidence of a light beam that is reflected by the sample and reaches the objectives. Thanks to its excellent technical features it is a equipment with a wide range of applications in industry, materials science, geology, etc.

#### Ergonomics

It is a very solid, large and stable equipment as well as with a comfortable and useful design. The head is 30° inclined and presents interpupillary distance adjustment that together with the diopter correction ring in one of the eyepieces allows an adequate observation. It is provided with a third tube for graphically recording images of the samples by using a digital photograph or video camera thus increasing considerably the equipment features. The light beam is directed by means of a sliding rod that sends 100% of the light to the eyepieces or to the third tube.

Sample focusing is performed through coarse and fine focusing knobs placed at both sides of the microscope in a lower position for a more comfortable handling. Moreover, the right knob is provided with a friction control ring and the left one with a focusing upper limit blocking lever to avoid the slide to accidentally hit the objectives.

#### Triple layer stage

The mechanical stage is very large (250x230 mm) and can be moved 154 mm along both axes. A special mechanism is used for moving the sample: the triple layer stage. It presents two movements in common with the double layer stages, but it is also provided with a third layer specialized in quick movements all along the stage, very useful for studying large samples.

#### Illumination

Zuzi metallurgical microscope model 101 is equipped with a Köhler illumination system that allows a better control over the light that reaches the sample. This illumination system includes a halogen lamp with intensity adjustment, a horizontally adjustable condenser and field and aperture diaphragms. Moreover, the lamp is provided with a centering mechanism to get a correct adjustment and thus assuring an optimum illumination and visualization of samples. The microscope also presents a filter holder with green, yellow, blue and matt glass filters to obtain a higher quality image.

#### Polarization

The microscope also allows the study of the optical properties of samples since it is equipped with an analyzer and a polarizer that can be rotated until polarizer and analyzer are orthogonally positioned. Both, analyzer and polarizer can be removed from the light path when they are not being used.

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