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**ASSEMBLY INSTRUCTIONS
FOR POLYPROPYLENE RACKS**

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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MAIN FEATURES

- Suitable for applications in aqueous media without loss of stability
- Resistant to autoclave sterilization processes
- Modular design allowing safe stacking of multiple racks
- Quick snap-fit system without the need for tools
- Numbering and layout facilitate sample identification
- Specific areas provided for the placement of labels or identification codes

ASSEMBLY PROCEDURE

1. Place the base on a flat surface with the word "BOTTOM" facing down.
2. Fold up one end plate and insert its lower tabs into the slots of the base. Make sure the imprinted numbers on the bottom tier face upward.
3. Position the bottom tier with the numbers facing upward and insert its tabs into the slots of the base and the end plate.
4. Fold up the second end plate and insert its tabs into the slots of the base and the bottom tier.
5. Insert the top tier into the upper slots of the end plates.
6. Fold up the side tier supports and insert their tabs into the slots of the base, the bottom tier, and the top tier.