BALANZA ELECTRÓNICA ELECTRONIC BALANCE BALANCE ÉLECTRONIQUE

Serie/Series/Séries 5161-LED

Ref. KBB012, KBB013

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita-Blue electronic balance, series 5161-LED. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and safety for its user.

The correct use of the equipment and its good quality will allow you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit shorts, fires, damages, etc. Please read the point of Tips, where we expose the safety notes.

<p>TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.</p>
--

Please bear in mind the following:

- This manual is inseparable from the Nahita-Blue electronic balance series 5161-LED, so it should be available for all the users of this equipment.
- You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the balance.
- To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- If you have any doubt about setting up, installation or functioning do not hesitate to contact your wholesaler.
- This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- Overhaul is not covered by the equipment warranty.
- Operations made by non-qualified staff will automatically void the warranty.
- Accessories (including their loss), are not covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- Manufacturer reserves the right to modify or improve the manual or equipment.

INDEX OF CONTENTS

1. SPECIFICATIONS	13
2. ASSEMBLY.....	14
3. TURN ON THE BALANCE.....	14
4. CALIBRATION.....	14
5. WEIGHTING.....	15
6. INTERFACE.....	16
7. TIPS.....	16
8. TROUBLESHOOTING.....	17
9. PACKING LIST.....	17

1. SPECIFICATIONS

Precision balances series 5161-LED work on high precision strain gauge load cell which implements high speed stabilization and high reliability.

Technical specifications

Code	KBB012	KBB013
Max. capacity	10kg	20kg
Tare range	10kg	20kg
Readability (g)	0.1g	0.1g
Repeatability (g)	+/-0.1g	+/-0.1g
Non-linearity (g)	+/-0.2g	+/-0.2g
Corner error (g)	+/-0.2g	+/-0.2g
Pan size (mm)	337x237mm	
Dimensions (LxWxH)	370x370x125mm	
Power supply	AC110V-220V Rechargeable battery 4V/4AH	

- Operating temperature range: 5-35°C
- Moisture: 50-85%

Features

- High precision load cell
- Rugged plastic housing
- Big size stainless steel pan
- High contrast easy to read LED display
- Built-in rechargeable battery
- Height adjustable feet
- Full capacity taring
- Fast response
- Overload protection
- Under balance weighing device
- Selectable measure units, kg, lb, g
- Weighting check, piece counting function
- RS232 interface

Function keys

ON/OFF, to turn on / off the balance

COU, for counting

TARE, for taring

UNIT, for unit conversion (g/lb/kg)

CAL, for calibration

MENU, for printing

2. ASSEMBLY

- Unpack the package, put the pan on the top of the loader.
- Put the balance on a steady flat surface away from vibrations, direct sunshine, air blows, strong magnetic disturbances and explosive, corrosive and flammable substances.
- Battery locker is at the rear of the balance, turn on the locker to turn on the balance. Please, turn off the locker during moving or transportation or if you are not going to use the balance for a long period of time.

3. TURN ON THE BALANCE

- Connect the balance to power supply, keep the switch in the ON(-) status.
- Press function key ON/OFF to turn on the balance, the balance will display following figures sequentially:

8.8.8.8.8.8.

Battery Voltage U=

Maximum capacity

- Finally the balance stays in a standby state showing 0g, 0.0g in the display

Notice: the lasting time of ----- will be decided by the stability of the load cell, thus, the balance must not be located in an unstable surface or exposed to wind blow. If O flashes in the display, it means the balance is not in a stable condition, calibration and counting operations are not allowed.

4. CALIBRATION

- The purpose for this operation is to calibrate the balance so as to achieve the best performance in case there is obvious tolerance error in weighting or the balance is under different gravity due to location in a different latitude.
- It is suggested to warm up the balance over half an hour before calibration.

- Remove all loads from the pan, press function key **TARE** to clear the readings to 0g, 0.0g.
- Press and hold function key **CAL** for 3 seconds, release function key **CAL** when “---CAL---“ is shown in the display, a figure of standard weight will flash in the display. Put a standard weight on the pan accordingly, standby state “-----“ will be shown in the display for a few seconds before the standard weight figure is shown in the display. Remove the weight, “-----“ standby state will remain for a few seconds before the balance enter stable weighing mode, showing 0g, 0.0g.
- It is recommended to make the operation twice to achieve the best calibration result.

5. WEIGHTING

When the balance is warmed up and calibrated, 0g, 0.0g is shown in the display indicating weighting mode. Put the object on the pan, when stable, the weight of the object is shown in the display.

Taring

When a container is put on the pan, its weight is shown in the display. Press function key **TARE**, 0g, 0.0g will be shown in the display, indicating the weight of the container is deducted. Put the object into the container, when stable, the figure shown in the display is the weight of the object.

Piece counting

Remove all loads from the pan, press function key **TARE** to clear the readings in the display, press function key **COU** to enter counting mode. Figure 10 (default sampling quantity) will be shown in the display, weighting unit changes from **g** to **pcs**. Put 10 samples on the pan, press function key **COU**, the balance will show “-----“ before it shows 10 pcs in the display. The balance enters counting mode.

Notice: In counting mode, the weight of the sample objects should be even, the weight of the individual sample should not be less than the readability of the balance.

Return to weighting mode

Press function key **COU** to return to weighting mode.

Unit conversion

Press function key **UNIT** to change weighing unit kg/lb/g, default weighing unit is set to g.

Overload

The weight of the object cannot exceed the rated maximum capacity of the balance, when exceeding, “-----“ will be shown in the display. Remove the object immediately from the pan so as to prevent damages to the balance.

- Always use original components and supplies. Other devices can be similar but they can damage the equipment
- If any liquid spills on the balance and comes in contact with the electric parts of the balance please do immediately disconnect it from power supply and send it to Nahita Technical Service as soon as possible for its checking and adjustment
- Never use the balance in a wedged location as for example a shelf.
- Never use sharp objects as pens, etc... to press buttons of the control panel; only use your fingers.
- Please, follow the instructions in this manual

8. TROUBLESHOOTING

Phenomenon	Possible Reason	Solution
Upper Line - - - -	Overload	Re-calibrate the balance
Under Line _ _ _ _	Overload or load cell broken	Re-calibrate the balance
Err-1	Too frequently turn on and off the balance	Turn off the balance, resume it after 3 seconds
Err-2	The balance is not stabilized	Wait for a few seconds for stabilization
	Low battery	Charge or replace battery

9. PACKING LIST

Description	QTTY
Balance	1 pcs
Pan	1 pcs
Power cord	1 pcs
Manual	1 pcs

NOTE: According to the in force legislation regarding “Non-automatic weighting instruments” in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 1/3rd/95), these balances must not be used for:

- Commercial transactions
- Calculating taxes, tariffs, rates, indemnities and other similar canons
- Judicial surveys
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories
- Determining the price or total amount in sale price or in pre-packaged preparations

Instructions on environmental protection

Nahita-Blue electronic balances series 5161-LED can operate with rechargeable batteries. Batteries contain heavy metals (zinc, lithium, mercury, etc.) that could contaminate water, soils, etc.

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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