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INDEX OF CONTENTS

Introduction and warranty	9	5. Operating procedure.....	13
1. Security instructions	10	6. Troubleshooting	13
2. Usage	10	7. Maintenance and cleaning	13
3. Inspection	11	8. Normative	14
4. Functioning	11	9. Technical specifications	14
4.1. Dashboard	11	10. Accessories	15
4.2. Display	12		

This is the user manual for the RSLAB-9 “roastery” and RSLAB-9 “disc” type stirrers. Please read this manual carefully before use. Instructions and procedures must be followed and the user must remain aware of the risks associated with the use of these devices.

If necessary, contact your dealer for help and advice on how to use the appliance correctly. To do this, the user must have the serial number of the device, describe the problems encountered, the procedures followed to try to solve them and the details of the contact person.

RSLAB stirrers are guaranteed for a period of 12 months from the invoice date against material and manufacturing defects under normal use conditions, as described in this manual. This warranty does not apply to the stirrer or any part damaged as a result of improper installation, improper connections, misuse, accident, or abnormal conditions of use.

Contact your dealer for any warranty claim. Returns are only possible with the agreement of the dealer.

1. SECURITY INSTRUCTIONS

!Attention!

Please read this manual carefully before using the appliance. Make sure that only qualified personnel use the appliance.

Protected electrical network

Make sure that the plug of the appliance is connected to the earth socket before using it.

Wear appropriate safety clothing for the risk category of the materials used, to avoid the risk of splashing and glass breakage.

Follow the safety instructions, guidelines and accident prevention regulations. To avoid risk of injury, do not touch live parts.

Place the appliance on a flat, stable, clean, dry and fireproof surface. Do not use the device in an explosive atmosphere, with hazardous materials or under water.

If the appliance does not work properly, reduce the speed.

Properly secure accessories and containers to prevent damage. Only use materials that do not react dangerously.

Use the standard accessories described in the Accessories section and follow the instructions for use of those accessories. Turn off the appliance when installing accessories and always check the appliance and accessories for damage before each use.

The device must only be opened by qualified personnel.

The voltage indicated on the appliance must correspond to that of the electrical network used. Do not cover the appliance and avoid shock to the appliance and its accessories.

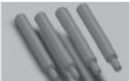
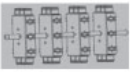
Keep the device out of the reach of a strong magnetic field.

2. USAGE

The apparatus has been designed to mix sticky materials in schools, laboratories or industry. It can be used with materials of different viscosities. The appliance is not suitable for domestic use or in an atmosphere that could be dangerous for the user or the appliance, as indicated in chapter 1.

The use of accessories not recommended by the manufacturer or the improper use of the appliance may endanger the user.

10. ACCESORIES

	Reference	Description
	LJJ007	60x centrifuge tubes Disc, 1,5 mL
	LJJ008	16x centrifuge tubes Disc, 15mL
	LJJ009	8x centrifuge tubes Disc, 50mL
	LJJ010	Adapter to assemble several discs
	LJJ001	Horizontal mounting for 48 x 1.5 mL centrifuge tubes
	LJJ002	Horizontal mounting for 24 x 15 mL centrifuge tubes
	LJJ003	Horizontal mounting for 24 x 50 mL centrifuge tubes
	LJJ004	Vertical mount of 32 centrifuge tubes of 1.5 mL
	LJJ005	Vertical mounting of 16 centrifuge tubes of 15 mL
	LJJ006	Vertical mounting of 16 centrifuge tubes of 50 mL

Before using any other cleaning or decontamination method, the user should consult the dealer to avoid damaging the device.

Wear suitable gloves to clean the appliance.

Do not use cleaning products to clean electronic components.

For any transport (for example, return for repair), the device must be cleaned, decontaminated and returned to its original packaging. In addition, suitable packaging must be used for transport.

If the appliance is not going to be used for a long period, unplug it and place it in a dry, clean place at a stable temperature.

8. NORMATIVE

The device complies with standard safety regulations:

EN61010-1, UL3101-1, CAN/CSA C22.2 (1010-1), EN61010-2-10

The device complies with EMC standards:

EN61326-1

Changes or modifications not expressly approved by authorities could void the user's authority to operate the equipment. This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These standards are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy. If not installed and used in accordance with the instructions in this manual, it may cause harmful interference to radio communications. Domestic use of this equipment may also cause harmful interference, in which case the user will be required to correct the interference at his own expense.

9. TECHNICAL SPECIFICATIONS

Tension	100-240 VCA
Frecuency	50/60 Hz
Power	40W
Rotation angle	0-90°
Motor type	CC Motor
Speed range	10-70 rpm
Speed display	LCD display
Timer	1-1199min
Timer display	LCD display
Operation type	Timed / Continuous
Outer dimensions	512x190x250mm
Weight	6,5 kg
Permissible ambient temperature	5-40°C
Permissible relative humidity	80%
Protection class DIN EN60529	IP21

3. INSPECTION

Carefully remove the device from its packaging and check that it has not been damaged during transport. If so, contact the dealer that supplied the appliance.

If the appliance is damaged, it must not be connected to the mains.

The RSLAB-9 “disk” and “roisserie” includes the following:

- Tube rotator, 1 unit
- Connection wire, 1 unit
- User manual, 1 unit

4. FUNCTIONING

4.1. Dashboard

Elements	Description
Power switch	On/Off
Angle control knob	Mando de control del ángulo de agitación.
LCD display	LCD screen showing real time and programmed values.
Speed/Timer light	According to the color of the light, the screen will show the speed or the duration.
Speed/Timer	When the light is green, the screen shows the speed. When the light is red, the screen shows the duration .
Power/Overload Protection Light Power/Overload	The light is green when the device is on. The light turns red when the overload protection system is activated. When the torque reaches its maximum value, the overload protection system is activated. At the same time it flashes and the appliance stops.
Mode change button	Toggles between displaying speed and time.
Speed control/timer knob	Select speed or duration by turning the knob. push on/off to activate/deactivate stirring .

4.2. Display

Parameter	Description
Set	Appears when values are programmed.
Err	Appears when an error occurs.
Display	When Set is displayed, the display shows the programmed value. When Set disappears, the display shows the real-time value.

5. OPERATING PROCEDURE

Place the shaker on a stable and firm surface and make sure that the voltage and power required for its use match the voltage and power of the mains. Make sure the power outlet is grounded. Plug in the appliance and turn it on. The appliance will start the self-diagnosis. After initialization is complete, the display shows the last speed value. Turn the speed and time selection knob to the desired speed.

- Press the value control button to start stirring. Press the button again to stop.
- Press the Speed/Timer mode change button and rotate the value control knob to set the desired time. Press the value control button to stop stirring. During operation, stirring can be stopped at any time by pressing the value control button. Pressing the button again will restart the agitation and run time will elapse. After the time has elapsed, the appliance stops automatically and emits an acoustic signal.

If all these operations are carried out normally, the appliance is ready for use. On the contrary, if any of these operations is not carried out normally, or if the appliance has been damaged during transport, contact your distributor for technical assistance..

Note:

The programmed values can be modified at any time.

6. TROUBLESHOOTING

The device does not turn on:

Check that the cable is correctly connected

Shaking stops suddenly:

If the overload protection light is red and the display shows "Er 03", the machine is overloaded. Press the value control button to stop the agitation function, reduce the load in the vessel and press the button again to resume agitation. If stirring does not resume, repeat the procedure, further reducing the vessel load.

7. MAINTENANCE AND CLEANING

Proper maintenance will keep your appliance in good condition and prolong its useful life. Make sure that no water or liquid enters the appliance when cleaning it. Unplug the appliance before cleaning it. Use the following cleaners:

- For dyes : Isopropyl alcohol
- For construction materials: Water + detergent / isopropyl alcohol
- For cosmetics: Water + detergent / Isopropyl alcohol
- For food : Water + detergent
- For fuels: Water + detergent