BALANZA ELECTRÓNICA SERIE 5162, 5161
ELECTRONICAL BALANCE SERIES 5162, 5161
BALANCE ÉLECTRONIQUE SÉRIES 5162, 5161

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Note: According to the in force legislation regarding “Non-automatic weighing instruments” in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 1/3rd/95), these balances must not be used for:

- Commercial transactions
- Calculating taxes, tariffs, rates, indemnities and other similar canons
- Judicial surveys
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories
- Determining the price or total amount in sale price or in pre-packaged preparations

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Instructions on environmental protection

Nahita-Blue electronic balances series 5161/5162 can operate with rechargeable batteries. Batteries contain heavy metals (zinc, lithium, mercury, etc.) that could contaminate water, soils, etc.

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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Spanish 2-8
 English..... 9-15
 French 16-22

INDEX OF CONTENTS

1. SPECIFICATIONS..... 10
 2. DESCRIPTION..... 11
 3. INSTALLATION..... 11
 4. USE OF THE BALANCE..... 13
 5. MAINTENANCE AND CLEANING 14

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita-Blue electronic balances, series 5161/5162. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years. The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- This manual is inseparable from the Nahita-Blue electronic balances 5161/5162, so it should be available for all the users of this equipment.
- You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the balance.
- To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler.
- This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- Overhaul is not covered by the equipment warranty.
- Operations made by non-qualified staff will automatically produce a loss of the warranty.
- Accessories (including their loss), are not covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage.
- In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- Manufacturer reserves the right to modify or improve the manual or equipment.

Overload

■ The weight of the sample cannot exceed the rated maximum capacity of the balance; when exceeded, the upper line "-----" will be displayed. Remove the sample immediately to avoid damages in the balance.

Energy saving mode

- Keep pressed COU key for 3 seconds to enter energy saving mode; press COU key and consecutively will be displayed:
 - CO-Off: display backlit is off
 - CO-1 : backlit of display will be off after 1 second of no use.
 - CO-5: backlit of display will be off after 5 seconds of no use.
- To exit energy saving mode, press CAL key successively until returning to weighing mode.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the balance it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- Please follow the processes of use of this manual.
- This manual should be available for all users of this equipment.
- Prevent the balance from sudden movements and knocks, as well as from direct sunlight or air flows. The balance is a precision instrument, you must handle it carefully.
- The balance is supplied with an adapter. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- Never unplug the adapter by pulling the wire, do it from the base.
- Never use the balance in a wedged location as for example a shelf.
- Never use sharp objects as pens, etc... to press the buttons of the control panel; only use your fingers.
- Never place on the pan an object heavier than the maximum capacity of the balance, the sensor could be damaged.
- Neither submerge the balance nor spill liquids on it.
- When you are not using the balance for a long period of time, lock the rechargeable battery.
- If any liquid comes into contact with the electric parts of the balance please do immediately disconnect it from the net and send it to the technical service, as soon as possible, for its checking and adjustment.
- Always use original components and supplies. Other devices can be similar but they can damage the equipment.

Cleaning

Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the balance and produce an early ageing of the equipment. Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

Note: If the balance has a series of discontinuous lines on the upper part of the screen, you must decide that your calibration has been lost due to transport or has a long time of disuse, to be able to eliminate them only requires a new calibration with the total weight.

Max. capacity	Calibration weight
300 g	200 g
500 g	500 g
1000 g	500 g
2000 g	1000 g
3000 g	2000 g
5000 g	2000 g

Before performing calibration, warm up the balance in advance.

- Keep pressed "TARE/CAL" key; The necessary weight value will flash.
- Place the calibration weight on the pan; Wait until the flashing weight value display changes to a steady state (e.g. 500.00 g).
- Remove the weight of the pan; "0.00" will be displayed and the balance will return to the weighing mode.

4. USE OF THE BALANCE

Weighing function

- Enter the weighing mode
- Place a container on the pan
- Once the weight on display has stabilized, press "TARE/CAL" key; value on display will change to 0.00.
- Place the sample to be weighed in the container.
- Once the value on display has stabilized, read the display.

Piece counting function

- Press "COU" key to enter in piece counting mode; "-COU-" will be displayed and then "10" (reference n° of pieces by default).
- Press "TARE/CAL" key to choose the desired reference number of pieces (10, 20, 50, 100, 150, 200, 250 or 500).
- Place the corresponding number of pieces on the pan and press COU key to save setting.
- Remove the pieces from the pan and the balance is ready to be used in piece counting mode.

Note: In counting mode, the weight of the pieces must be even and the weight of an individual piece should not be less than the readability of the balance.

- To return to weighing mode, press "COU" key.

Weighing unit selection

- Balance can perform weighing in different measuring units.
- In weighing mode press repeatedly "UNIT" key to consecutively display all available weighing units and select the desired one.
- The weighing unit by default is g.

1. SPECIFICATIONS

Precision balances series 5161 and 5162 work on high precision strain gauge load cell which implements high speed stabilization and high reliability.

SERIES 5161


Code	KBB025	KBB004	KBB005	KBB006	KBB007
Max. capacity (g)	500	1000	2000	3000	5000
Tare range (g)	0-500	0-1000	0-2000	0-3000	0-5000
Readability	0.1g				
Pan size (mm)	120*155mm				
Dimensions (LxAxH)	280x180x80mm				
Power supply	AC 110-120V/220-240V				

SERIES 5162

Code	KBC001	KBC002	KBC003	KBC004
Max. capacity (g)	300	500	1000	2000
Tare range (g)	0-300	0-500	0-1000	0-2000
Readability	0.01g			
Pan size (mm)	120*155mm			
Dimensions (LxAxH)	280x180x80mm			
Power supply	AC 110-120V/220-240V			

- High precision load cell
- Full plastic housing
- Stainless steel pan
- Easy to read LCD with backlit
- Overload display indicator
- Low or charged battery display indicator
- Mains adapter supplied as standard
- With built-in rechargeable battery
- Lock/Unlock battery switch
- Height adjustable feet
- Full capacity taring
- Simple 4 key menú for easy operation
- Automatic calibration
- Selectable measuring units: g, kg, ct, T, TAR, dr, PKT, GN, TMR, gsm, tIJ, mo, dwt, oz, lb, tIT, ozt, tIH, %.
- Piece counting function

2. DESCRIPTION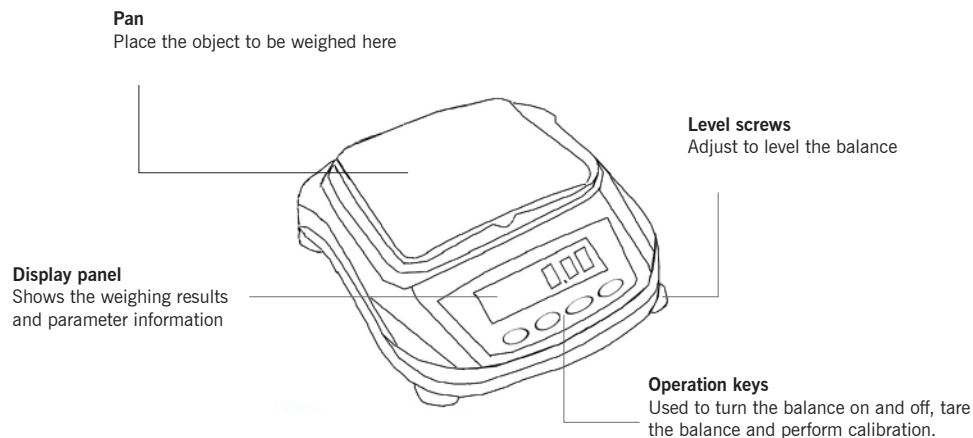

Operation keys function

Key	In weighing mode	
	Press once and release	Keep pressed for a while
ON/OFF	Turn the balance on and off	-
COU	Select the piece counting function	(Always after entering the calibration menu) Linear calibration
TARE/CAL	Tare function (set weighing value to 0)	Calibration function
UNIT	Select the weighing unit	-

3. INSTALLATION

Choosing the installation place

- The measuring performance of the balance is greatly influenced by the environment where it is installed.
 - Observe the following points to ensure a safe and accurate weighing.
 - Do not use the balance anywhere exposed to explosive, combustible or corrosive gases. This could cause fire or other kind of serious problems.
 - Use the correct power supply and voltage with the balance. Using an incorrect power supply or voltage with the balance will lead to fire or trouble with the balance.
- Note also that if the power supply or voltage is unstable or if the power supply capacity is insufficient, it will not be possible to obtain satisfactory performance from the balance.
- Avoid locations where the balance will be exposed to any of the following situations:
 - Air flow from an air conditioner, ventilator, door or window.
 - Extreme temperature changes

- Vibration from surroundings or nearby equipments.
- Direct sunlight.
- Dust, electromagnetic waves or a magnetic field.

■ Install the balance on a strong and stable flat table or floor. Placing the balance in an unstable site could lead to troubles with the balance or injury to users. When selecting the installation site, take into account the combined weight of the balance and the item to be weighed.

Unpacking and delivery inspection

■ Check that all of the items indicated below are included in the package, and that nothing has been damaged:

- Balance
- Pan
- Instruction manual
- Adapter


■ In case the balance or any component presents any damage due to transport tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Turning the power on

- Insert the plug of the power cord into the DC IN connector at the back of the balance.
- Unlock the rechargeable battery by the switch at the bottom of the balance.

Note: if you are not using the balance for a long period of time lock the battery to preserve it in good conditions.

- Connect the power cord to the power outlet.
- Press "ON"; The display will automatically light on and go through the changes indicated below: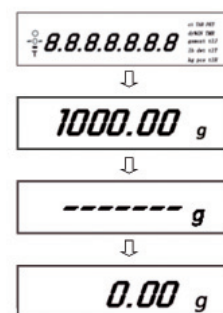

Warming up

- Before performing calibration on the balance or measuring its accuracy it is necessary to ensure that the balance is in a stable state.
- For balance stabilization it is important that environmental temperature is stable.
- Leave the balance on and in weighing mode for at least half an hour before performing calibration; this is called "warming up".

Calibration

- Perform balance calibration each time after moving it. Weights are required for external calibration balances.
- Necessary calibration weight depending on maximum capacity of the balance