DECLARACION DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DECLARATION DE CONFORMITE CE

El fabricante | The manufacturer | Le fournisseur:  
**AUXILAB, S.L.**

Declara que el equipo | Declare that the equipment | Declara que l'appareil:

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Directive for electromagnetic compatibility (EMC)

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Electrical equipment for measurement, control and laboratory use. EMC requirements.

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Safety requirements for electrical equipments for measurement, control and laboratory use. Part 1: General requirements.

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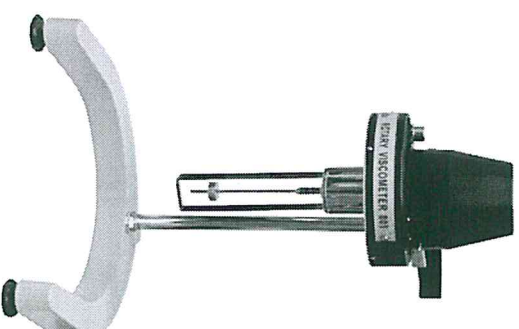
This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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- ◆ Never try to repair the viscosimeter by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the viscosimeter (burns, hurts...) or damages in nearby equipments.

◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

#### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the viscosimeter and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.
- ◆ For cleaning spindles, use neutral products with pH not higher than 6.

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**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

#### **INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita viscosimeter 801. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita viscosimeter 801, so it should be available for all the users of this equipment.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the viscosimeter.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Accessories (including their loss), are not covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

base of current 220 V, 50 Hz  $\pm$  10%

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220V 50Hz  $\pm$  10%.

#### Setting up

- ◆ Remove the protecting cap that covers the thread of the spindle holder (1.6) and screw the selected spindle by turning it clockwise (to unscrew, turn the spindle in anticlockwise direction).

**NOTE: Hold the spindle holder (1.6) with one hand and with the other carefully screw the spindle.**



- ◆ Prepare the fluid to be measured in a beaker or a cylindrical container of at least 70 mm diameter and control its temperature precisely.
- ◆ Slowly, raise the recipient with the fluid until the spindle is submerged up to the circular mark.
- ◆ Release the indicator braking lever (1.5) and switch the motor on by pressing power on/off (1.2); turn the speed adjustment knob (1.4) until the desired speed is upwards matching with the red mark. After several turns, 20 or 30seconds, indicator reading tends to be stable. Readings can also be made at specific times.
- ◆ If the rotating speed is fast (30 or 60 r.p.m.) press the indicator braking lever (1.5) to cause the reading to settle.
- ◆ When the indicated value is too high or too low, change spindle or speed; it is better to perform measurements at a speed range from 30-90 rpm.

#### Selection of range, coefficient, spindle and speed

- ◆ First of all, the approximate range of viscosity of the fluid must be estimated. Then, select the appropriate speed and spindle according to table of ranges (Chart 1).
- ◆ For instance, if the estimated viscosity is 3000 cP, one of the following combinations must be selected: spindle n° 2 at a speed of 6 r.p.m or spindle n° 3 at a speed of 30 r.p.m.
- ◆ When viscosity is difficult to estimate, a high viscosity must be supposed. Try the spindle from larger to smaller and speed from lower to higher, until finding the most adequate combination for the viscosity of the fluid.

CHART 1. According to speed and spindle

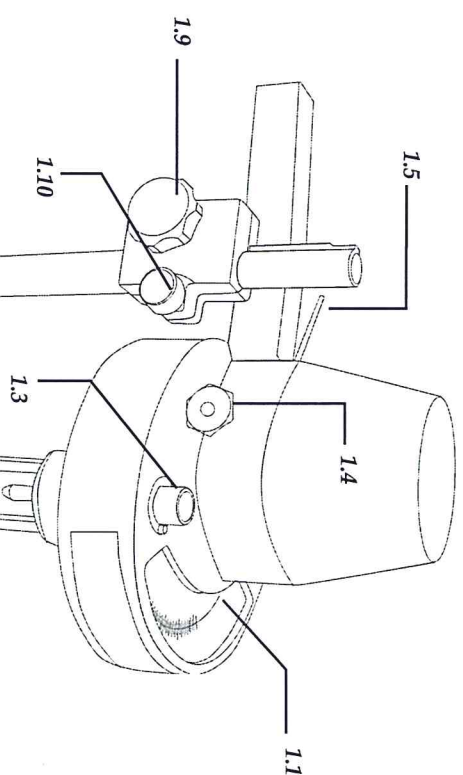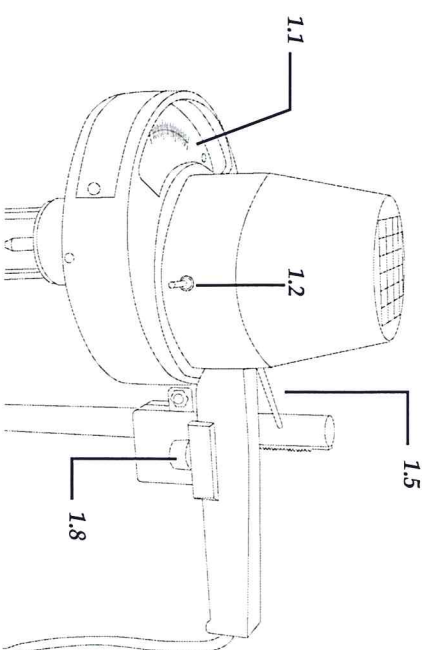Spindle	60	30	12	6
1	100	200	500	1000
2	500	1000	2500	5000
3	2000	4000	10000	20000
4	10000	20000	50000	100000

### 3. TECHNICAL SPECIFICATIONS

Measuring range	10-100000 mPa.s (cP)
Precision	1%
Spindles	N° 1, 2, 3 and 4
Rotor speed	6, 12, 30 and 60 r.p.m.
Tolerance	$\pm$ 5% (on Newtonian fluids)
Power	220 V, 50-60 Hz