SET DE CUBOS DE METAL
METAL CUBE SET
SET DE CUBES EN MÉTAL

Nahita

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EXPERIMENT 2**DETERMINE WHETHER AN OBJECT WILL FLOAT OR SINK IN WATER**

Extra material required:

- Beaker with water

Procedure:

1. Take the beaker with water
2. Insert the different metal cubes into the water
3. Observe the results.

Table:

Material/Objeto	Flota/Hunde

Result:

Materials that sink in water have a higher density than water, and materials that float in water have a lower density than water.

Applications:

Density can be used to calculate the weight of a piece of machinery or a part of a bridge or construction before it is built. An engineer only needs to know the volume and density of the material it is made of. The engineer can pre-calculate the load that each part of the structure will bear. This will determine whether the design is sturdy enough.

Density can be used for other types of pure metals and alloys, as in the classic case of Archimedes where he had to determine whether the king of Sicily had a crown made of pure gold or alloy.

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INDEX OF LANGUAGES

Spanish 2-4
 English 5-7
 French 8-10

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Description

Cubes for density investigation available in different materials: brass, lead, iron, copper, aluminum, and zinc. These cubes are used for learning about density, mass, volume, specific gravity, and flotation phenomenon. They are also used for verifying the Archimedes principle.

Density:

All matter has mass and volume. These are physical properties of materials and vary depending on the material. The amount of matter contained in a material is called mass, and the amount of space occupied is called volume. Density is the measure of mass per unit of volume.

$$\text{Density} = (\text{mass of a substance} / \text{volume of a substance})$$

Where D is the density of an object and V is the volume of the object.
 In SI units, density is expressed in kg/m³.
 For example, the density of water is 1000kg/m³

Material	Densidad del material (g/ml)
Copper	8.9
Brass	8
Steel	7.6
Aluminum	2.7
Acrylic	1.16-1.19
Oak	0.60-0.90
Nylon	1.13
Pine	0.35-0.60
Polar	0.35-0.05
PVC	1.39-1.42
Polypropylene	0.85-0.95
Tree of life	1.28-1.37

EXPERIMENT 1

DETERMINE THE DENSITY OF A GIVEN OBJECT USING A SPRING BALANCE

Extra material required:

- Spring balance or dynamometer.
- Beaker with water
- Cable

We can measure the density of an object by the weighing method. This is one of the most accurate methods for determining it.

1. Suspend the sample using a cable from the spring balance and record the weight.
2. Lift the beaker with water around the sample, completely submerging the sample, and re-weigh the sample.
3. The difference between the weights is the weight of water displaced. From that value, and knowing the density of water (1gm/ml)

$$\text{Density} = (\text{Weight of sample in air}) * (\text{density of water}) / (\text{Weight of sample in air} - \text{Weight of sample in water})$$

Table:

Material	Weight in air (P1, grams)	Weight in water (P2, grams)	P1-P2 (grams)