## Vasos comunicantes

Referencia: QLB004

### Descripción

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Code: QLB004

### Description

The apparatus comprises of four glass tubes, each of different cross-sectional size and shape, connected together at the bottom through a common horizontal glass tube, sealed to a manifold and mounted on stand. It demonstrates that the liquid level in communicating vessels remains same irrespective of the shapes and sizes of the vessels.

### Communicating vessels

Set of containers containing a homogeneous fluid. When the liquid settles it balances out to the same level in all of the containers. If additional liquid is added to the vessel, the liquid will again find a new equal level in all the connected vessels. The process of equal level water in different vessels is a part of Stevin's Law.

### Stevin's Law

It states that the pressure at a point in a liquid in static equilibrium depends on the depth at that point.

$$P = P_o + dgh$$

Where

$P_o$  is the external pressure on the surface of the liquid

$h$  is the depth

$g$  is the gravity acceleration

$d$  is the liquid density

All the vessels contain equal level of water because gravity and pressure are constant in each vessel.

### Experiment

To demonstrate that the liquid level in communicating vessels remains same irrespective of the shapes and sizes of the vessels.

### **Other components required**

- Beaker
- Funnel

### **Procedure**

1. Fill colored water in the beaker.
2. Put the funnel on the mouth of any one of the vessels and pour the water in the vessel from the beaker.
3. Fill the Apparatus at least half level of the vessels.
4. Remove the funnel from the vessels and observe that the water takes the shapes of the vessels with the same levels.

### **Conclusions**

From experiments, we conclude that water does not a definite shape and it take the shape of the different shapes of the vessel in which is water is filled. The water level remains the same when vessels of the different shapes and sizes are connected to each other.

### **Precautions**

- Performs the experiments carefully since the apparatus is fragile in nature.
- Always fill the colored liquid in the vessels.